

References:

1. Luczynska CM, Arruda LK, Platts-Mills TA, Miller JD, Lopez M, Chapman MD. A two-site monoclonal antibody ELISA for the quantification of the major Dermatophagoides spp. allergens, Der p I and Der f I. J Immunol Methods 1989; 118(2):227-235.
2. Custovic A, Taggart SC, Francis HC, Chapman MD, Woodcock A. Exposure to house dust mite allergens and the clinical activity of asthma. J Allergy Clin Immunol 1996; 98(1):64-72.
3. Platts-Mills TA, Vervloet D, Thomas WR, Aalberse RC, Chapman MD. Indoor allergens and asthma: report of the Third International Workshop. J Allergy Clin Immunol 1997; 100(6 Pt 1):S2-24.
4. Sawyer G, Kemp T, Shaw R et al. Biologic pollution in infant bedding in New Zealand: high allergen exposure during a vulnerable period. J Allergy Clin Immunol 1998; 102(5):765-770.
5. Arbes SJ, Jr., Cohn RD, Yin M et al. House dust mite allergen in US beds: results from the First National Survey of Lead and Allergens in Housing. J Allergy Clin Immunol 2003; 111(2):408-414.
6. Gruchalla RS, Pongracic J, Plaut M et al. Inner City Asthma Study: relationships among sensitivity, allergen exposure, and asthma morbidity. J Allergy Clin Immunol 2005; 115(3):478-485.
7. Zock JP, Heinrich J, Jarvis D et al. Distribution and determinants of house dust mite allergens in Europe: the European Community Respiratory Health Survey II. J Allergy Clin Immunol 2006; 118(3):682-690.
8. Jarvis D, Zock JP, Heinrich J et al. Cat and dust mite allergen levels, specific IgG and IgG(4), and respiratory symptoms in adults. J Allergy Clin Immunol 2007; 119:697-704.
9. Earle CD, King EM, Tsay A et al. High-throughput fluorescent multiplex array for indoor allergen exposure assessment. J Allergy Clin Immunol 2007; 119(2):428-433.



www.inbio.com

Indoor Biotechnologies, Inc.
700 Harris Street,
Charlottesville, VA 22903
United States

Indoor Biotechnologies Ltd.
Vision Court
Caxton Place
Cardiff, Wales CF23 8HA
United Kingdom

Indoor Biotechnologies,
India Private Limited
Bangalore Bioinnovation Centre,
BioTech Park, Electronic City Phase 1
Bangalore-560100, India

Tel: (434) 984-2304
Fax: (434) 984-2709
Email: mail@inbio.com

Tel: +44 (0) 29 2167 4640
Email: Info@indoorbiotech.co.uk

Tel: +91-9901722009
Email: info@inbioindia.com




Der p 1 ELISA kit (10B9/5H8)

Product Code: EL-DP1A

Lot Number: XXXXX

Sample Curve:

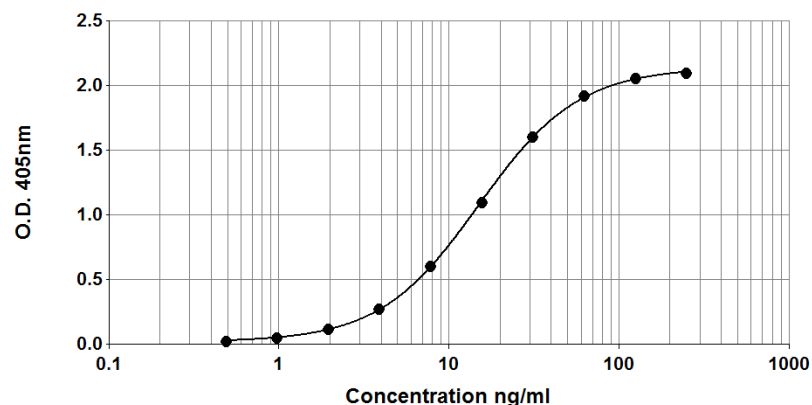

Content:

Vial 1 (red top) 100 μ l
Monoclonal antibody 10B9

Vial 2 (white top) 400 μ l
Universal Allergen Standard
Concentration: 2500ng/ml Der p 1

Vial 3 (brown) 100 μ l
Biotinylated monoclonal antibody 5H8
Dilute: 1:1000 for use

Storage: The ELISA kit should be stored at 4°C

For research and commercial use in vitro: not for human in vivo or therapeutic use.

An InBio™ product

Certificate of Analysis

Monoclonal Antibody: 10B9 (clone 10B9 F6 A12)
Immunogen: Der p 1
Isotype: Mouse IgG1
Specificity: Binds to an epitope on dust mite *Dermatophagoides pteronyssinus* allergen, Der p 1.
Purification: Produced in ascites and purified by ammonium sulphate precipitation and affinity chromatography using Protein A. Single heavy and light chain bands on SDS-PAGE.
Concentration: 1.0mg/ml in phosphate buffered saline, pH 7.4. Based on A280 for IgG (1.42= 1mg/ml), 0.22µm filtered, preservative free.

Lot Number: xxxxx

Monoclonal Antibody: 5H8 (clone 5H8 C12 D9)
Immunogen: Der p 1
Isotype: Mouse IgG2A
Specificity: Binds to an epitope on dust mite *Dermatophagoides pteronyssinus* allergen, Der p 1.
Purification: Produced in cell culture and purified by affinity chromatography using Protein A. Single heavy and light chain bands on SDS-PAGE.
Biotinylation: Biotinylated and titrated for use in ELISA at 1/1000 dilution. Prepared in 1% BSA/50% glycerol/PBS, pH 7.4, 0.22µm filtered, preservative free.

Lot Number: xxxxx

Allergen Standard: Universal Allergen Standard
Composition: A formulation of eight purified natural allergens prepared in 1% BSA/50% glycerol/PBS, pH 7.4

Lot Number: xxxxx

Concentration:

Universal Allergen Standard	Protein Measurement	Concentration (ng/ml)
Der p 1	Amino-acid analysis	2500
Der f 1	Amino-acid analysis	2500
Der p 2	Amino-acid analysis	1000
Fel d 1	Amino-acid analysis	1000
Can f 1	Amino-acid analysis	2500
Rat n 1	Amino-acid analysis	1000
Mus m 1	Amino-acid analysis	250
Bla g 2	Amino-acid analysis	2500

ELISA protocol for Der p 1.

- Coat polystyrene microtiter plates (NUNC Maxisorp Cert. NUNC catalog # 439454) with 100µl mAb 10B9 at 10µl/10ml, i.e. 1/1000 dilution of stock, in 50mM carbonate-bicarbonate buffer, pH 9.6, incubate overnight at 4°C.
- Wash wells 3x with PBS-0.05% Tween 20, pH 7.4 (PBS-T). Incubate for 30 min. at room temperature with 100 µl/well of 1% BSA, PBS-T. Wash 3x with PBS-T.
- Use doubling dilutions of the Universal Allergen Standard to make a control curve ranging from 250 - 0.5ng/ml Der p 1: Pipette 20µl UAS into 180µl 1% BSA, PBS-T into wells A1 and B1 on the ELISA plate. Mix well and transfer 100µl across the plate into 100µl 1% BSA, PBS-T diluent to make 10 serial doubling dilutions. Wells A11, B11 and A12, B12 should contain only 1% BSA, PBS-T as blanks.
- Add 100µl of diluted allergen samples and incubate for 1 hour at room temperature. House dust extracts for Der p 1 analysis are routinely diluted two-fold from 1/10-1/80. Other sample types, like air filter extracts and allergen extracts, may require different dilutions.
- Wash wells 3x with PBS-T and add 100µl diluted biotinylated anti-Der p 1 mAb 5H8. The antibody solution contains 50% glycerol and should be diluted 1/1000 in 1%BSA, PBS-T. Incubate for 1 hour at room temperature.
- Wash wells 3x with PBS-T and add 100µl diluted Streptavidin-Peroxidase (Sigma S5512, 0.25mg reconstituted in 1ml distilled water). The reconstituted Streptavidin should be diluted 1/1000 in 1% BSA, PBS-T. Incubate for 30 minutes at room temperature.
- Wash wells 3x and develop the assays by adding 100µl 1mM ABTS in 70mM citrate phosphate buffer, pH 4.2 and 1/1000 dilution of H₂O₂. Read the plate when the absorbance at 405nm reaches 2.0-2.4.

Notes:

Universal Allergen Standard is recommended for immunoassay calibration purposes only. Not recommended for in-vitro antibody measurements, T cell studies, immunization purposes, or other uses.

Buffer recipes, storage conditions and a list of frequently asked questions can be found under "Protocols" on our web site: www.inbio.com.

For research and commercial use in vitro: not for human in vivo or therapeutic use.